


CLEANING AND PUTTING AWAY THE OLYMPUS CH30 MICROSCOPE

1. Make sure all oil is removed from the oil-immersion lens with lens paper.
2. Use a clean (non-oily) piece of lens paper to clean off the 10X lens.
3. Put the 10X objective in place.
4. Wrap cord around the base of the microscope.
5. Put the microscope into the cabinet with the handle facing outward, so it can be pulled out easily by the next user.

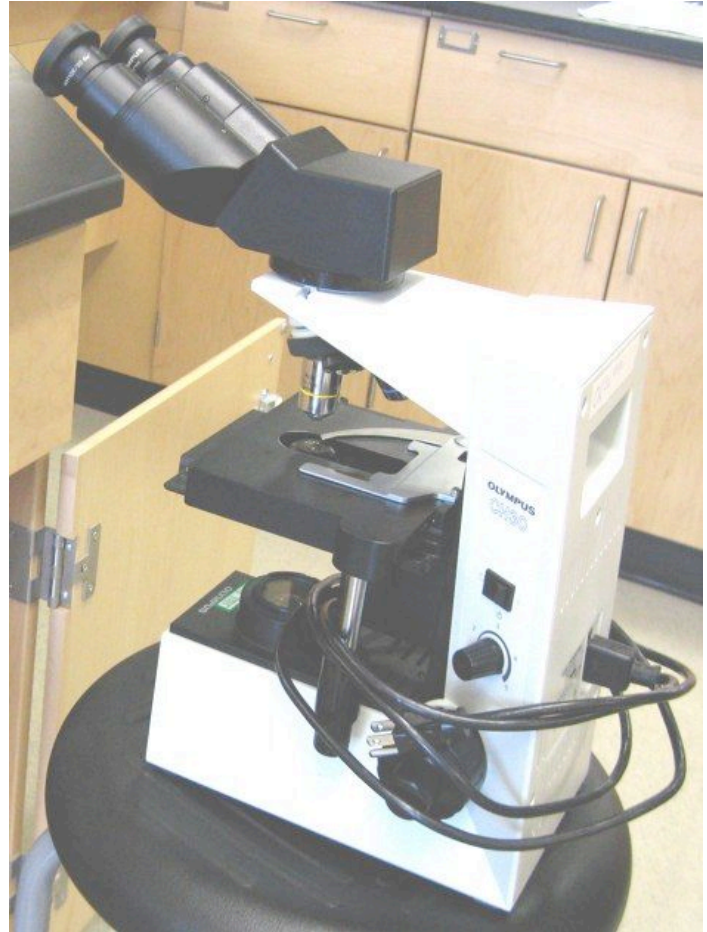

USING THE PHASE-CONTRAST MICROSCOPE

(for viewing wet mounts only)

PRELIMINARIES:

1. Have the 40X objective lens in place.
2. Place slide on stage such that it is secured along three sides. (Don't put slide under the clip.)
3. Please do not pull this unit out! This is what makes the phase scope work to bring out the necessary contrast.

FOCUSING:

4. With the coarse adjustment, bring the slide up very close to the objective lens while looking from the side.
5. While looking through the scope, use the fine adjustment to pull the image into focus.

